
## Data Logger/Acquisition Software

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Seite 2–5

**GB** Instruction Manual

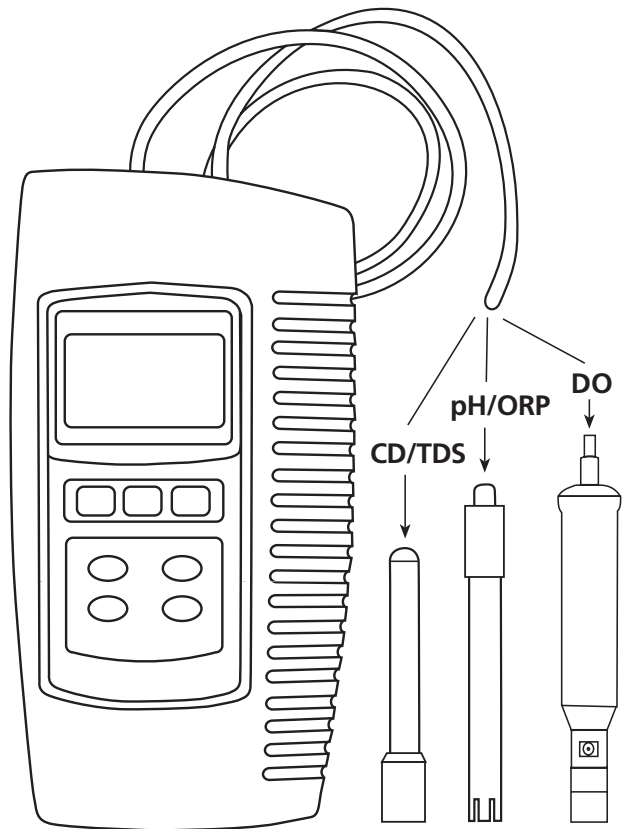Page 6–8

---

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---

|    |                                      |   |
|----|--------------------------------------|---|
| 1. | <b>Allgemeine Beschreibung</b> ..... | 3 |
| 2. | <b>Softwareinstallation</b> .....    | 3 |
| 3. | <b>SW-E802.xls</b> .....             | 3 |
| 4. | <b>Meter_Test</b> .....              | 5 |

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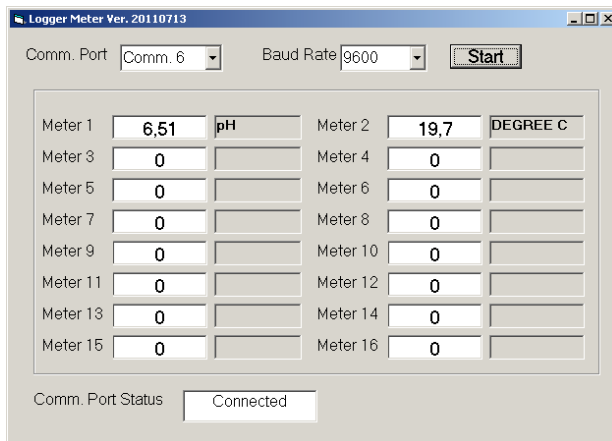
## 1. General Description

The SW-E802 application software is designed to transfer and log the data generated by the AL15 meter in an Excel Spread Sheet. Also an on-line display of the instrument's data on a PC is possible.

An RS232 cable is required to transfer the data. Should the computer not have an RS232 interface, the instrument also can be connected via a USB cable. A virtual COM port can be established using the drivers supplied on the CD.

Part No.: 724500 (RS232 cable)

Part No.: 724510 (USB cable with virtual COM port)

## 2. Software Installation

If an older version of the software has previously been installed, please delete it before installing the new version. To install the software, insert the CD into the CD ROM disk drive of the computer. After the automatic start-up, please follow the instructions of the installation wizard.

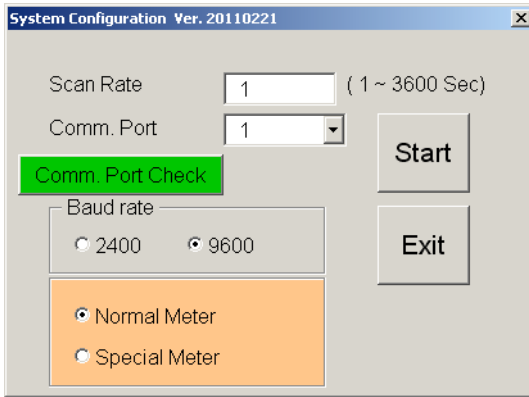
On successful installation, a new programme SW-E802 will appear the "Start/All programs" directory.

## 3. SW-E802.xls

On execution of the SW-E802.xls, the following Excel-Sheet will appear on the screen.

|   | A              | B         | C                 | D         | E        | F         | G        | H         | I        | J         | K        |
|---|----------------|-----------|-------------------|-----------|----------|-----------|----------|-----------|----------|-----------|----------|
|   | Config / Start | Stop Date | Manual Input Time | Ch1 Value | Ch1 Unit | Ch2 Value | Ch2 Unit | Ch3 Value | Ch3 Unit | Ch4 Value | Ch4 Unit |
| 1 |                |           |                   |           |          |           |          |           |          |           |          |
| 2 |                |           |                   |           |          |           |          |           |          |           |          |
| 3 |                |           |                   |           |          |           |          |           |          |           |          |
| 4 |                |           |                   |           |          |           |          |           |          |           |          |
| 5 |                |           |                   |           |          |           |          |           |          |           |          |
| 6 |                |           |                   |           |          |           |          |           |          |           |          |

Click the 'Config/Start' button to access the below screen where all the required settings can be selected. Confirm the settings with the 'Start' button. The data logging will now begin.


The following settings may be selected:

- Select the scan rate.
- Choose the COM port used to connect the hand held meter. 'Comm. Port Check' enables the user to select the required COM interface.
- Set the baud rate (recommended: 9600).
- Select the meter (Normal meter).
- Select the 'Exit' button to close the setting window without saving the changes.
- Click the 'Start' button to save the settings and start the data logging.

|   | A              | B          | C                 | D         | E        | F         | G        | H         | I        | J         | K        | L   |
|---|----------------|------------|-------------------|-----------|----------|-----------|----------|-----------|----------|-----------|----------|-----|
|   | Config / Start | Stop Date  | Manual Input Time | Ch1_Value | Ch1_Unit | Ch2_Value | Ch2_Unit | Ch3_Value | Ch3_Unit | Ch4_Value | Ch4_Unit | Ch5 |
| 1 |                |            |                   |           |          |           |          |           |          |           |          |     |
| 2 | 1              | 05.12.2011 | 08:54:53 5,58     | pH        | 19,7     | DEGREE C  |          |           |          |           |          |     |
| 3 | 2              | 05.12.2011 | 08:54:58 5,58     | pH        | 19,7     | DEGREE C  |          |           |          |           |          |     |
| 4 | 3              | 05.12.2011 | 08:55:03 5,58     | pH        | 19,7     | DEGREE C  |          |           |          |           |          |     |
| 5 |                |            |                   |           |          |           |          |           |          |           |          |     |
| 6 |                |            |                   |           |          |           |          |           |          |           |          |     |
| 7 |                |            |                   |           |          |           |          |           |          |           |          |     |

The data will continue to be logged until the 'Stop' button is selected.

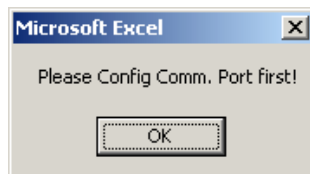The 'Manual Input' button enables the user to store single readings to the Excel-sheet. Transmission is only possible if the settings have been correctly selected. Otherwise the following message appears.


The 'Clear' button can be selected to delete all data logged on the Excel-Sheet. The following warning window will appear. Select 'Ja' to confirm deletion of data or 'Nein' to cancel.

#### 4. Meter\_Test

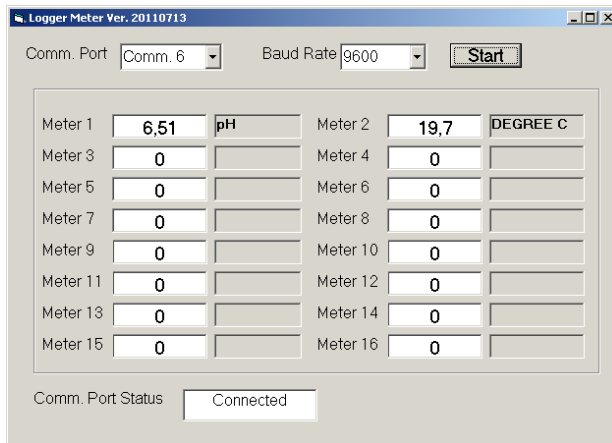On starting Meter\_Test, the following operating window is displayed. Once the correct COM port for connection and the baud rate (recommended: 9600) have been selected, the reading of the hand held meter can be displayed by clicking the 'Start' button.

---

---

---

**Tintometer GmbH, Division Aqualytic®**

Schleefstraße 8-12  
44287 Dortmund  
Tel.: (+49) (0)2 31 / 9 45 10-755  
Fax: (+49) (0)2 31 / 9 45 10-750  
sales@aqualytic.de  
www.aqualytic.de  
Germany

Technical changes without notice  
Printed in Germany 12/11