Photometer AL250

Chlor HR • Chlorine HR • Chlore HR
Cloro HR • Cloro HR

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Seite 3–14

(GB) Instruction Manual

Page 15–26

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Page 27–38

(IT) Istruzioni d'uso

Pagina 39–50

(ES) Instrucciones

Página 51–61

**CE-Konformitätserklärung / Declaration of CE-Conformity
Déclaration de conformité CE / Dichiarazione di conformità CE /
CE-Declaración de conformidad**

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- **General notes** 16
 - Guidelines for photometric measurements 16
 - Method notes 17

- **Functional description** 18
 - Operation 18
 - Measuring the absorbance 19

- **Methods** 20
 - Chlorine HR, with tablet 20

- **Menu options** 22
 - Menu selections 22
 - Recall of stored data 22
 - Setting date and time 22

- **Calibration Mode** 23
 - User calibration 23
 - Recommended calibration values 23
 - Factory calibration reset 24


- **Technical data** 25
 - Operating messages 25
 - Error codes 25

 **CAUTION** 

The accuracy of the instrument is only valid if the instrument is used in an environment with controlled electromagnetic disturbances according to DIN 61326. Wireless devices, e.g. wireless phones, must not be used near the instrument.


Guidelines for photometric measurements

1. Vials, caps and stirring rods should be cleaned thoroughly after each analysis to prevent interferences. Even minor reagent residues can cause errors in the test result.
2. The outside of the vial must be clean and dry before starting the analysis. Clean the outside of the vials with a towel. Fingerprints or other marks will be removed.
3. Zero calibration and test must be carried out with the same vial as there may be slight differences in optical performance between vials.
4. The vials must be positioned in the sample chamber for zeroing and test with the ^l mark on the vial aligned with the Δ mark on the instrument.
5. Always perform zeroing and test with the vial cap tightly closed.
Place the cover on the adapter.
6. Bubbles on the inside wall of the vial lead to incorrect measurements. To prevent this, remove the bubbles by swirling the vial before performing the test.
7. Avoid spillage of water into the sample chamber because this can lead to incorrect test results.
8. Contamination of the lens in the sample chamber can result in errors. Check at regular intervals and – if necessary – clean the light entry surfaces of the sample chamber using a moist cloth or cotton buds.
9. Large temperature differences between the instrument and the environment can lead to errors – e.g. due to the formation of condensation in the area of the lens or on the vial.
10. To avoid errors caused by stray light do not use the instrument in bright sunlight.
11. Always add the reagent tablets to the water sample straight from the foil without touching them with the fingers.
12. The reagents must be added in the correct sequence.

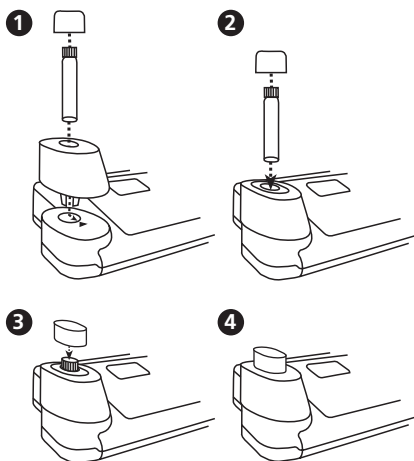Correct position of the vial (Ø 16 mm):**Method notes**

- Prior to measurement ensure that the sample is suitable for analysis (no major interferences) and does not require any preparation i.e. pH adjustment, filtration etc.
- Reagents are designed for use in chemical analysis only and should be kept well out of the reach of children.
- Ensure proper disposal of reagent solutions.
- Material Safety Data Sheets: available on request
(Internet: www.aqualytic.de)

GB Functional description

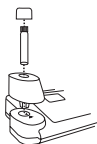CLHr

METHOD

METHOD

0.0.0

METHOD

RESULT

Operation

Fix the adapter for 16 mm vials on the sample chamber.

Switch the unit on using the [ON/OFF] key.

The display shows the following:

Select the required test using the [MODE] key:
CLHr → Abs1 → CLHr →..... (Scroll)

The display shows the following:

Fill a clean vial (16 mm Ø) with 8 ml of the water sample, close tightly with the cap. Place the vial in the adapter making sure that the Δ marks are aligned. Place the cover on the adapter.

Press the [ZERO/TEST] key.

The "Method" symbol flashes for approx. 8 seconds.

The display shows the following:

After zero calibration is completed, remove the vial from the sample chamber. The characteristic coloration appears after the addition of the reagents.

Close the vial tightly with the cap. Place the vial in the adapter making sure that the Δ marks are aligned. Place the cover on the adapter.

Press the [ZERO/TEST] key.

The "Method" symbol flashes for approx. 3 seconds.

The result appears in the display.

The result is saved automatically.

Repeating the test:

Press the [ZERO/TEST] key again.

New zero calibration:

Press the [MODE] key until the desired "Method" symbol appears in the display again.

Display backlight

Press the [!] key to turn the display backlight on or off. The backlight is switched off automatically during the measurement.

Recall of stored data


If the instrument is switched on, press the [!] key for more than 4 seconds to access the recall menu.

Measuring the absorbance

Abs

This Mode function allows measurement of the absorbance at the specified wavelength. The result appears in the display in mAbs. 1000 mAbs = 1 A (Absorbance unit)

Abs1: $\lambda = 470 \text{ nm}$

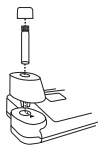0.0.0

Chlorine HR with Tablet 5 – 200 mg/l

Fill a clean vial (16 mm Ø) with **8 ml of the water sample** and perform zero calibration (see "Operation").

Remove the vial from the sample chamber.

Add **one CHLORINE HR (KI) tablet** straight from the foil to the water sample and crush the tablet using a clean stirring rod.

Add **one ACIDIFYING GP tablet** straight from the foil to the same water sample and crush the tablet using a clean stirring rod.

Close the vial tightly with the cap and swirl several times until the tablets are dissolved.

Place the vial in the adapter making sure that the Δ marks are aligned. Place the cover on the adapter.

Press the [ZERO/TEST] key.

The method symbol flashes for approx. 3 seconds.

The result is shown in the display in mg/l Chlorine.

CLHr

RESULT

Tolerance: ± 5 mg/l

Notes:

1. Oxidising agents interfere as they react in the same way as Chlorine.

Mode

On
Off

!

Mode

Menu selections

Press the [MODE] key and **hold**.

Switch the unit on using the [ON/OFF] key.

Allow the 3 decimal points to be displayed before releasing the [MODE] key.

The [!] key allows for selection of the following menu points:

- ▲ recall stored data
- ▲ ▾ setting the date and time
- ▾ user calibration

The selected menu is indicated by an arrow in the display.

Confirm the selection with the [MODE] key.

▲ Recall of stored data

The photometer shows the last 16 data sets in the following format (automatically proceeds every 3 seconds until result is displayed):

Number	n xx (xx: 16...1)
Year	YYYY (i.e. 2008)
Date	mm.dd (monthmonth.dayday)
Time	hh:mm (hourhour:minute)
Test	Method
Result	x,xx

The [ZERO/TEST] key repeats the current data set.

The [MODE] key scrolls through all stored data sets.

Quit the menu by pressing [!] key.

Zero
Test

Mode

!

2 3 Setting date and time (24-hour-format)

After confirming the selection with the [MODE] key the value to be edited will be shown for 2 sec.

The setting starts with the year (YYYY) followed by the actual value to be edited. Same applies for month (mm), day (dd), hour (hh) and minutes (mm). Set the minutes first in steps of 10, press the [!] key to continue setting the minutes in steps of 1.

Increase the value by pressing the [MODE] key.

Decrease the value by pressing [ZERO/TEST] key.

Proceed to the next value to be edited by pressing [!] key.

After setting the minutes and pressing the [!] key the display will show "IS SET" and instrument returns to the measurement mode.

ATTENTION: If the battery is removed for more than one minute the date and time menu starts automatically when the photometer is switched on the next time.

Mode

SET

DATE

YYYY

(2 sec.)

Mode

Zero
Test

!

4 User calibration

Note:

user calibration (Display in calibration mode)
factory calibration (Display in calibration mode)

CAL

CAL

CAL

CLHr

After confirming the selection with the [MODE] key the instrument will show CAL/CLHr.

Scroll through methods using the [MODE] key.

Fill a clean vial with 8 ml of the standard, close tightly with the cap and place the vial in the adapter making sure that the X marks are aligned. Place the cover on the adapter.

METHOD

0.0.0

CAL

Press the [ZERO/TEST] key.

The method symbol flashes for approx. 8 seconds.

The display shows the following in alternating mode:

Perform calibration with a standard of known concentration (see "Operation").

Press the [ZERO/TEST] key.

METHOD

RESULT

CAL

The method symbol flashes for approx. 3 seconds.

The result is shown in the display, alternating with CAL.

If the reading corresponds with the value of the calibration standard (within the specified tolerance), exit calibration mode by pressing the [ON/OFF] key.

Changing the displayed value:

Pressing the [MODE] key once increases the displayed value by 1 digit.

Pressing the [ZERO/TEST] key once decreases the displayed value by 1 digit.

CAL

RESULT + x

On
Off

Press the corresponding key until the reading equals the value of the calibration standard.

By pressing the [ON/OFF] key, the new correction factor is calculated and stored in the user calibration software.

:

Confirmation of calibration (3 seconds).

Recommended calibration values

Chlorine HR:

between 70 and 150 mg/l

Factory calibration reset

Resetting the user calibration to the original factory calibration will reset all methods and ranges.

A user calibrated method is indicated by an arrow while the test result is displayed.

To reset the calibration press both the [MODE] and [ZERO/TEST] key and **hold**.

Switch the unit on using the [ON/OFF] key.

Release the [MODE] and [ZERO/TEST] keys after approx. 1 second.

The following messages will appear in turn on the display:

SEL

CAL

The factory setting is active.
(SEL stands for Select)

or:

SEL

CAL

Calibration has been set by the user.

(If the user calibration is to be retained, switch the unit off using the [ON/OFF] key).

Mode

SEL

CAL

Calibration is reset to the factory setting by pressing the [MODE] key.

The following messages will appear in turn on the display:

On
Off

Switch the unit off using the [ON/OFF] key.

Technical Data

Light source:	LED, Filter ($\lambda = 470 \text{ nm}$)
Battery:	9 V-block battery (Life approx. 600 tests, without display light)
Auto-OFF:	Automatic switch off 5 minutes after last keypress
Ambient conditions:	5–40°C 30–90% rel. humidity (non-condensing)
CE:	Certificate for Declaration of CE-Conformity at www.aqualytic.de


To ensure maximum accuracy of test results, always use the reagent systems supplied by the instrument manufacturer.

Operating messages

Hi	Measuring range exceeded or excessive turbidity.
Lo	Result below the lowest limit of the measuring range.
	Replace 9 V battery, no further tests possible.
btLo	Battery capacity is too low for the display backlight; measurement is still possible.

Error codes

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E 01	Light absorption too great. Reasons: e.g. dirty optics.
E 10 / E 11	Calibration factor "out of range"
E 20 / E 21	Too much light reaching the detector.
E 22	Battery capacity was too low during measurement. Change battery.
E 70	CL HR: Factory calibration incorrect / erase
E 71	CL HR: User calibration incorrect / erase

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Tintometer GmbH, Division Aqualytic®